

## Portable LCD Digital Microscope LCD.62.G6



### Main parameters:

- 1: Pixel: HD 3 .6MP CCD
- 2: Display: 4.3 inches HD OLED display.
- 4: Magnification: 1-600X continuous magnification system.
- 5: Object distance: 15mm-- infinity
- 6: With built-in lithium battery, Continuous working for 6 hours .
- 7: Languages: English, Spanish, Russian, Korean,


Japanese, Thai, Hebrew, Portuguese, German, French, Italy, Turkey, Czech, Poland, Traditional Chinese, simplified Chinese.



The High-capacity lithium battery needs about 4 hours to be full charged. When it's charging, the display will show

🔌, when it's full charged, the display will show 🔋

(Important Note :if shipped by Air, Lithium battery is prohibited , we will change it as AC adapter Power Supply)

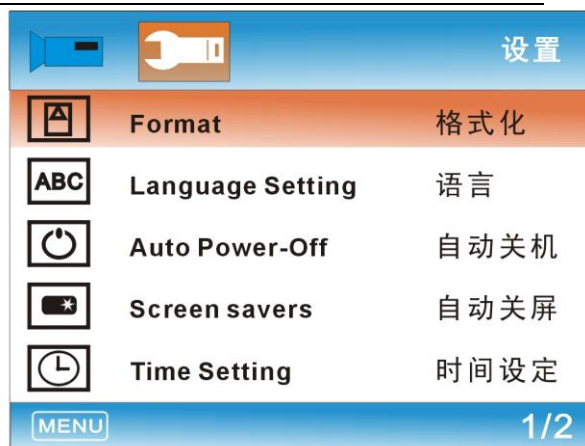

## : **Power On/Off**

Press it once to turn on, press once again to turn off.

## : **Function setting key:**

1. Resolution: 1080P, 720P, VGA
2. Date Tag: Displayed /Not displayed
3. Motion Detection: ON/OFF (When it's on it will start to record a video automatically , as soon as detecting something's moving .)
4. Cyclic Record : off /3 minutes /5 minutes /10 minutes.  
In the memory card is full, it will delete the first video and save the new video automatically.
5. Exposure value: +2.0 (0) -2.0
6. Format: Yse / NO
7. Language: 16 kinds of languages


8. Automatic shutdown: off, 1 minutes, 3 minutes,
9. Automatic turn off the LCD : off, 1,3,5 minutes
10. Factory Reset: Yes / No
11. Light source's frequency: 50HZ, 60HZ
12. Time setting: year, month, day, hours, minutes, seconds
13. Image rotating on / off
14. Light: on / off
15. Software version: V1.10

### Mode button:

1. Recording Video mode:
2. Taking Photos mode: (5M, 2M, 1.3M, VGA) press "OK" button for taking photos
3. Playback mode: delete, lock / unlock




Left button, right button, OK button (photo / video, open key)

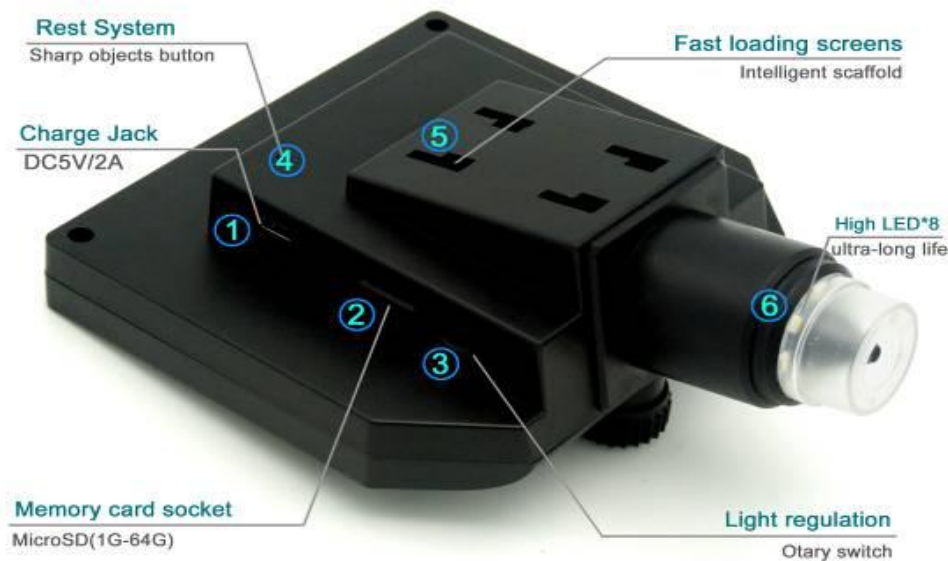

- 1: Power DC interface (5P MINI).
- 2: Memory card's socket (Micro SD), supporting 1-64GB.
- 3: Brightness Adjust button
- 4: REST button, restore the system.(If the machine can not work, please press the button with sharp objects to restart it.)
- 5: Fast loading 4 holes' bayonet, push the bracket's 4 claws into the bayonet,hearing "ka" once,it says the assemble is finished .
- 6: High brightness 8 LEDs, available for up to 100 thousand hours.

### **Intelligent universal support**




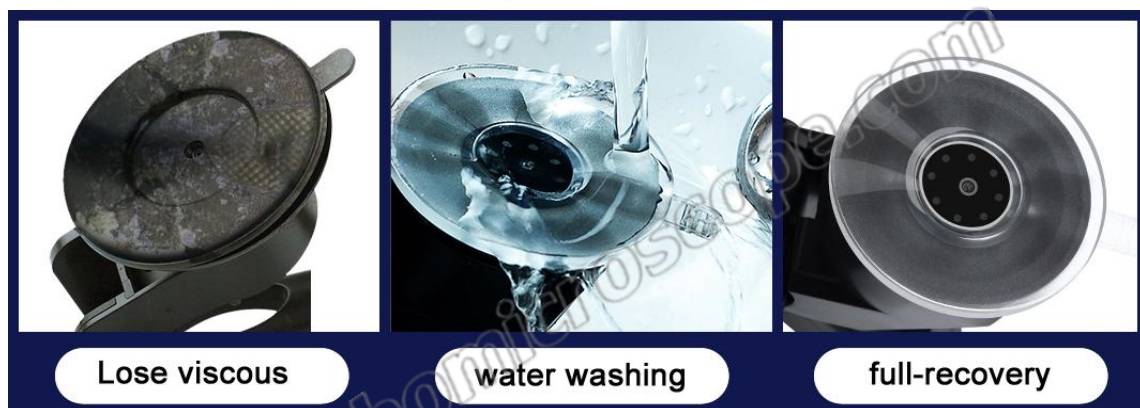

The intelligent universal support is super light and powerful, can be adsorbed on the surface of glass/ceramic/marble/plastic board and others.

The silicone sucker will not leave marks on any surface. after long-term using, it can be cleaned by water when there's dust on the surface.

The closest distance of microscope and the goal is 15mm, the magnifying time is the maximum , the magnifying time will be smaller as the distance increasing .

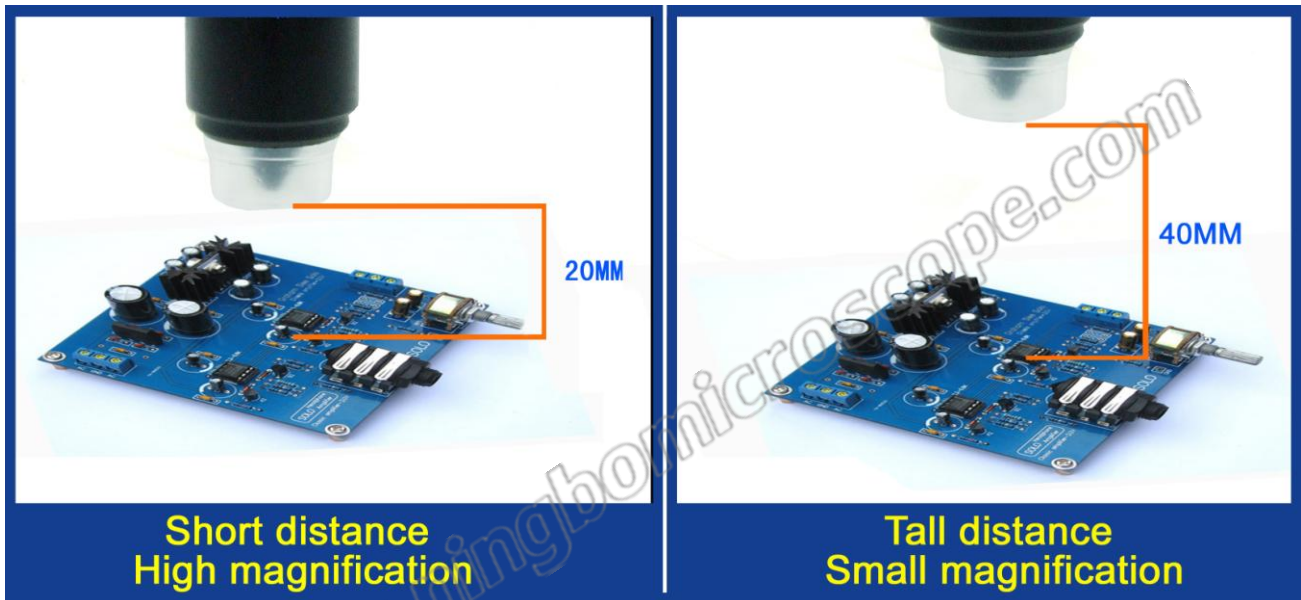

## Accessories icon:



Microscope, power supply, data line, multi functional support, user's manual.